

Lab Instructions


AOAC EXTRACTION METHOD

- 1** •Check Expiry Date
•Check Desiccant Pouch:
Blue = OK
Pink = Do not use test




2

Place indicated volume of buffer in vial A and vial B.

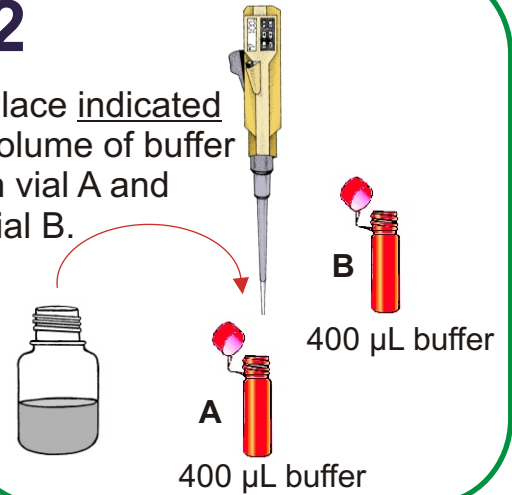


3

Add 100 µL of shellfish extract to 1.5 mL of water and mix. To mix, cap vial and shake thoroughly.

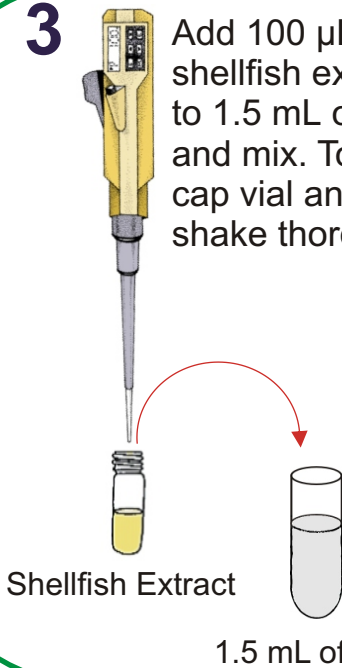

4


Take 100 µL of the shellfish extract/water mixture. Add to vial A and mix.

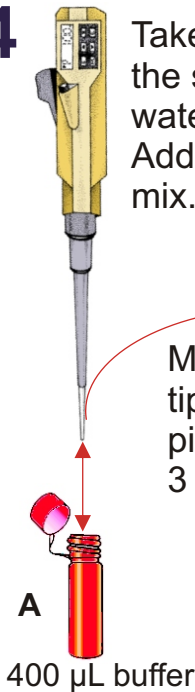

5

Take 100 µL from vial A. Add to vial B and mix.


Mix by inserting pipette tip into liquid and pipetting at least 3 times.

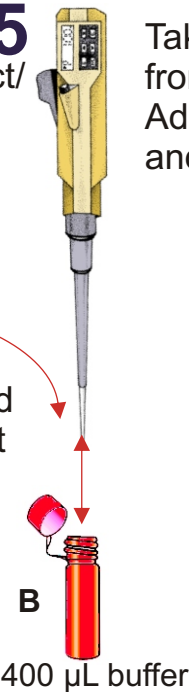


6

Place 100 µL of mixture from vial B into the sample well "S".

DO NOT REUSE PIPETTE TIPS

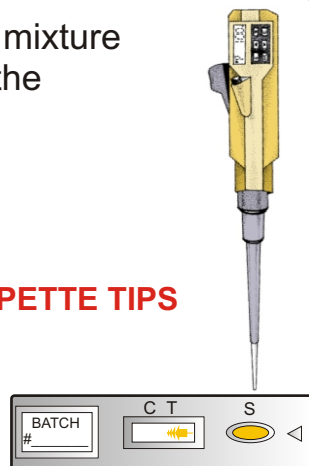

7

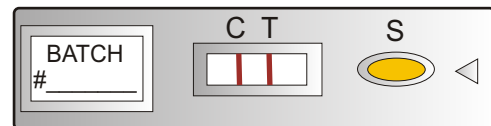

Read your results after 45-75 minutes.

Turn page for C & T Line Interpretation.