



# Lab Instructions METHANOL EXTRACTION METHOD


- 1** •Check Expiry Date  
•Check Desiccant Pouch:  
Blue = OK  
Pink = Do not use test

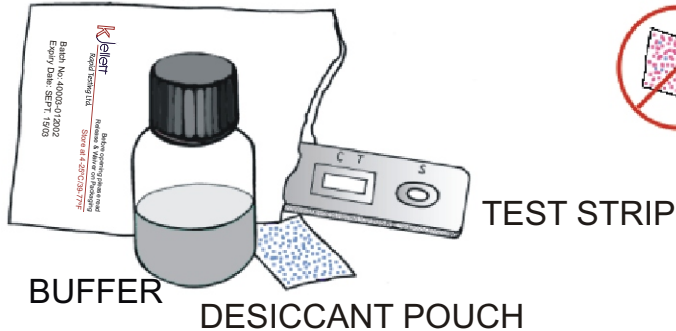


- 2** Place 900  $\mu$ L of buffer in a vial.

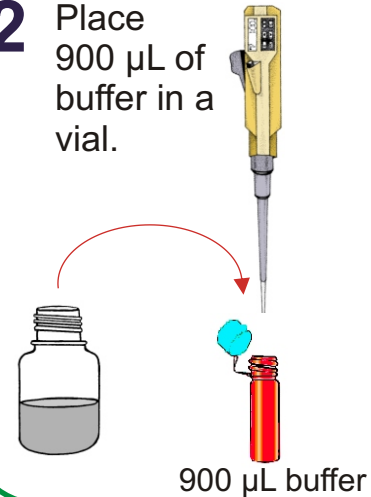

- 3** Add 100  $\mu$ L of shellfish extract to 1.5 mL of water and mix. To mix, cap vial and shake thoroughly.

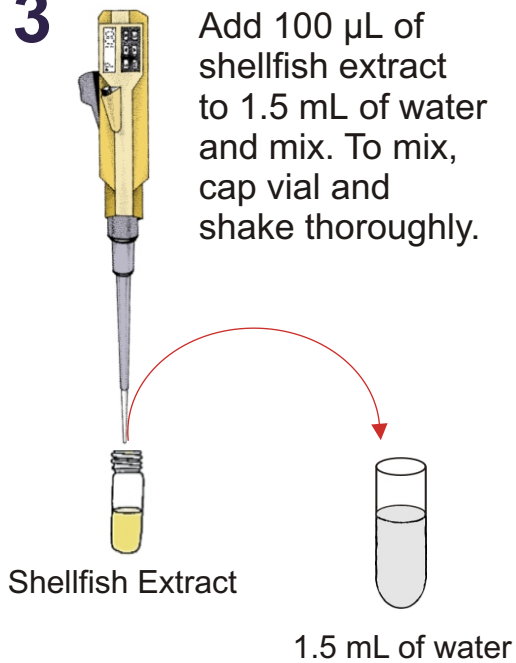


- 4** Take 100  $\mu$ L of the shellfish extract/water mixture to the vial in step 2 and mix.

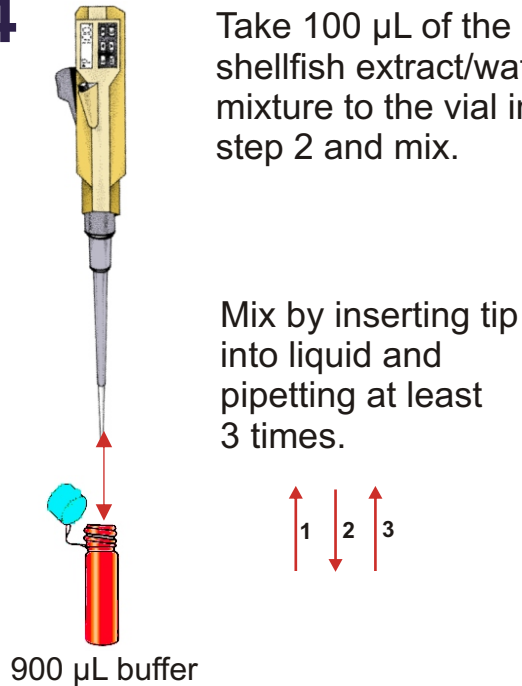


- 5** Place 100  $\mu$ L of mixture from the vial in step 4 into the sample well "S".

**DO NOT REUSE PIPETTE TIPS**

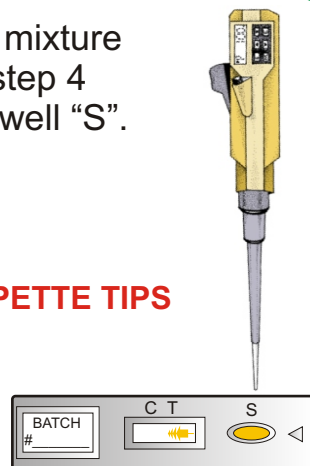

**6**

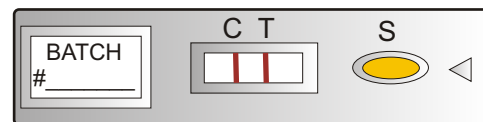

Read your results after 45-75 minutes.

Turn page for C & T Line Interpretation.