


# Field Instructions

# PHYTO EXTRACTION METHOD

- 1**
- Check Expiry Date
  - Check Desiccant Pouch:  
Blue = OK  
Pink = Do not use test

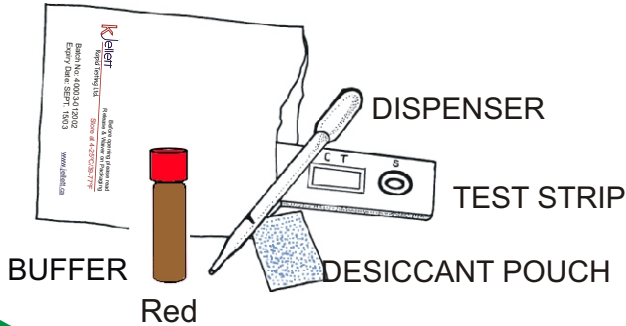

- 2** Before using, hold buffer vial as shown, 1 inch above table.




Tap 3 times on a table to get the liquid from the lid to the vial.


1"

Table

- 3** Add 100  $\mu$ L of phyto extract (to black line of dispenser) to the vial in step 2.

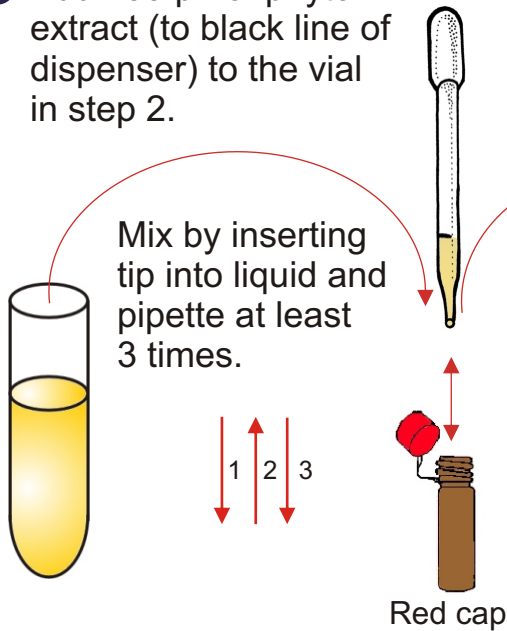

- 4** Place 100  $\mu$ L of the mixture into sample well "S".

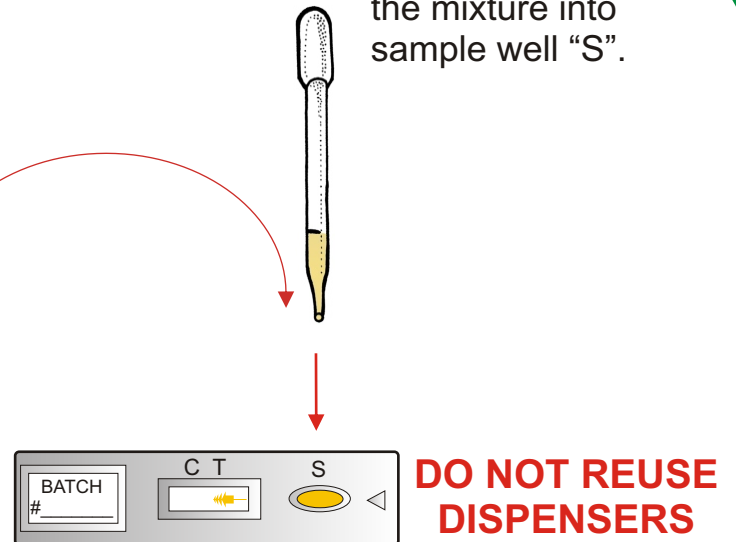

**5**

Read your results between 45-75 minutes.  
Turn this page for C & T Line Interpretation.

